

Sharing Data with Users Offsite

1. Setup Completed by Users

1. Login to globus.org using CLS credentials. This needs to be completed prior to sharing data as this will create a user account on Globus.

Log in to use Globus Web App

Use your existing organizational login

e.g., university, national lab, facility, project

Canadian Light Source

Didn't find your organization? Then use [Globus ID to sign in](#). ([What's this?](#))

Continue

Or

Sign in with Google

Sign in with ORCID iD

2. Copying Data to DTN Node (Beamline Staff)

3. Sharing Folder with Users (Beamline Staff)


4. User Data Access

1. From endpoints, select shared with you. This will show any collections shared with user. Select "REIXS User Data Access".

The screenshot shows the Globus Endpoints interface. The top navigation bar includes a search bar for endpoints and a link to create a personal endpoint. Below the navigation bar, there are filter tabs: 'Recently Used', 'In Use', 'Shareable by You', 'Shared with You' (which is highlighted with a blue box), and 'Administered by You'. A dropdown menu for the 'Shared with You' filter shows 'Filter shared with you'. The main content area is a table with columns: ENDPOINT, STRICT, STATUS, ROLE, and SHARED. The table contains one entry: 'REIXS User Data Access' with a status of 'ready'. Below the table name, it says 'Shared Endpoint on Canadian Light Source DTN'. The left sidebar contains navigation icons for File Manager, Bookmarks, Activity, Endpoints (selected), Groups, Console, Account, Logout, and Help.

ENDPOINT	STRICT	STATUS	ROLE	SHARED
REIXS User Data Access Shared Endpoint on Canadian Light Source DTN		ready		

2. Select "Open in File Manager".

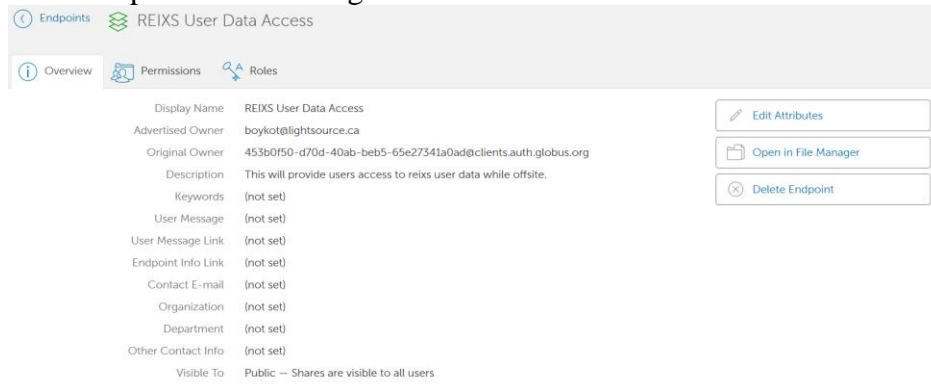

Endpoints REIXS User Data Access

Overview Permissions Roles

Display Name REIXS User Data Access [Edit Attributes](#)

Advertised Owner boykot@lightsources.ca [Open in File Manager](#)

Original Owner 453b0f50-d70d-40ab-beb5-65e27341a0ad@clients.auth.globus.org [Delete Endpoint](#)

Description This will provide users access to reixs user data while offsite.

Keywords (not set)

User Message (not set)

User Message Link (not set)

Endpoint Info Link (not set)

Contact E-mail (not set)


Organization (not set)

Department (not set)

Other Contact Info (not set)

Visible To Public — Shares are visible to all users

3. Navigate to your desired file, select and click "Download". You will be asked re-authenticate your account.

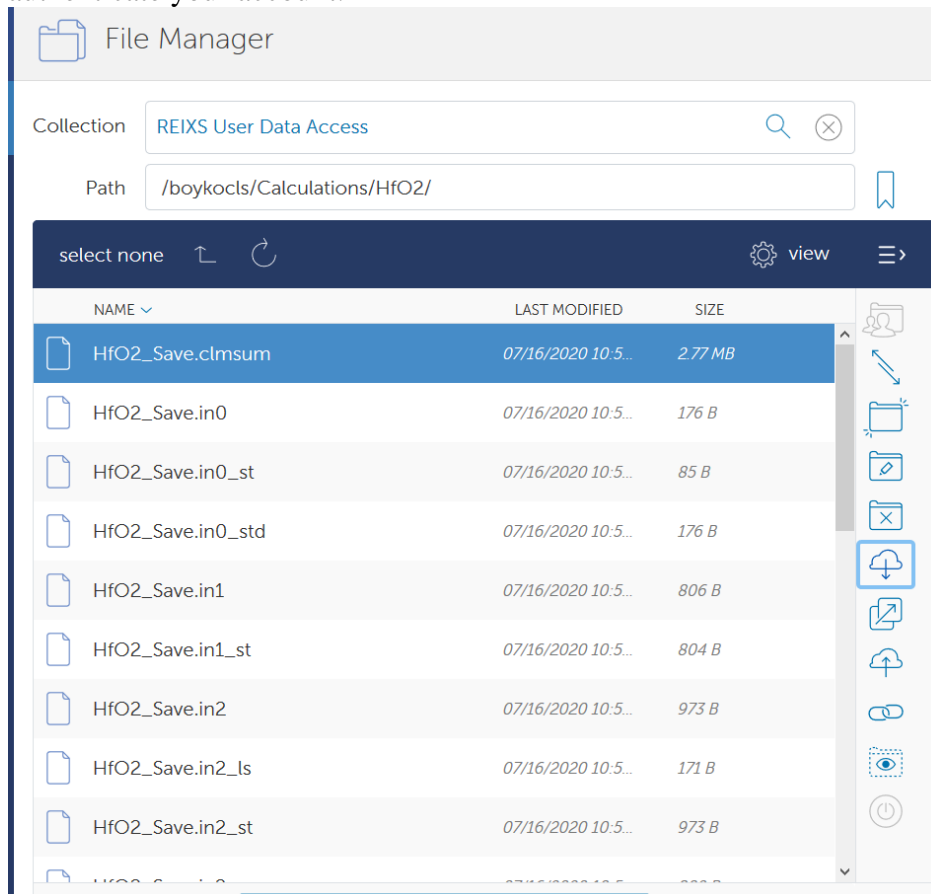

File Manager

Collection REIXS User Data Access

Path /boykocls/Calculations/HfO2/

select none ↶ ↷ view ≡

NAME	LAST MODIFIED	SIZE
HfO2_Save.clmsum	07/16/2020 10:5...	2.77 MB
HfO2_Save.in0	07/16/2020 10:5...	176 B
HfO2_Save.in0_st	07/16/2020 10:5...	85 B
HfO2_Save.in0_std	07/16/2020 10:5...	176 B
HfO2_Save.in1	07/16/2020 10:5...	806 B
HfO2_Save.in1_st	07/16/2020 10:5...	804 B
HfO2_Save.in2	07/16/2020 10:5...	973 B
HfO2_Save.in2_ls	07/16/2020 10:5...	171 B
HfO2_Save.in2_st	07/16/2020 10:5...	973 B